



<b>Composition</b>	Acetic Acid 3% in Purified Water.
<b>Use</b>	An acidic solution, which decreases the pH of the surrounding environment in laboratory applications.
<b>Warnings</b>	Do not irrigate the brain, meninges or perforated eardrum. The solution is not isotonic. The solution is haemolytic.
<b>Adverse Effects</b>	A transient stinging sensation may occur if spilt on skin.
<b>Directions for Use</b>	Use as necessary to acidify surrounding media.
<b>Safety Directions</b>	Avoid contact with eyes and skin. Avoid breathing vapour.
<b>First Aid</b>	For advice, contact a Poisons Information Centre (phone Australia 131126) or a doctor. If swallowed, do NOT induce vomiting.  If in eyes, hold eyelids apart and flush the eye continuously with running water. Continue flushing until advised to stop by the Poisons Information Centre or doctor, or for at least 15 minutes.
<b>Storage</b>	Store below 25°C
<b>Poison Schedule</b>	Not scheduled.
<b>Presentation</b>	

Code	Product	Size	Carton
ACE01214F	ACETIC ACID 3% v/v	100mL	20

**Contact Point**

Drug Information Pharmacist  
ORION Laboratories Pty Ltd ABN 56 009 293 136  
25-29 Delawney Street, Balcatta, Western Australia 6021 AUSTRALIA  
**Telephone (all hours): +618 9441 7800**  
FREE PHONE: 1800 805 546 FREE FAX: 1800 004 110  
EMAIL: customerservice@orion.net.au; WEBSITE: www.orion.net.au  
ORION® is a registered trademark of Orion Laboratories Pty Ltd

Prepared by: Anna McLean	Date: July 2011
Checked by: Robert Kimpton	Date: July 2011

**Revision Information**

Date	Document Name	Superseded Document	Revision Information
June 2006	ACE01214_2 June 2006	ACE01214_1 July 2003	General review; change of address
June 2009	ACE01214_03 June 2009	ACE01214_2 June 2006	General review
November 2009	ACE01214_04 November 2009	ACE01214_03 June 2009	<i>First Aid</i> : inclusion of phone number
July 2011	ACE01214_05 July 2011	ACE01214_04 November 2009	General review